


3.3 Upgrade Installation

A Upgrade Installation is when you run a installer that is newer than the version of IonWizard that is currently installed.



If you want to install an old (earlier) version that what is currently installed you must remove the current (newer) version and then do a normal installation of the older version.


To start the installer **double-click** on the MSI file from the Windows explorer (My Computer)

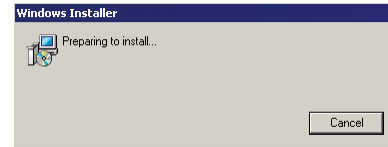

If you are upgrading a version of IonWizard that was installed in c:\ionoptix\current you will see a dialog that explains you are upgrading a "pre-msi" version.




After upgrading a pre-msi installation you should NOT try to run IonWizard from c:\ionoptix !

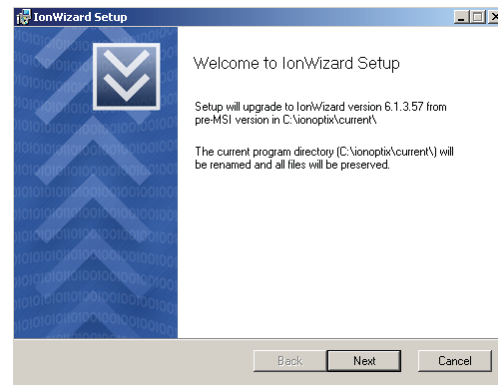

If you are upgrading a version of IonWizard that was installed in c:\ionoptix\current you will see the software license agreement

Check the box to confirm that you accept the license terms then **click** the Next button to continue.

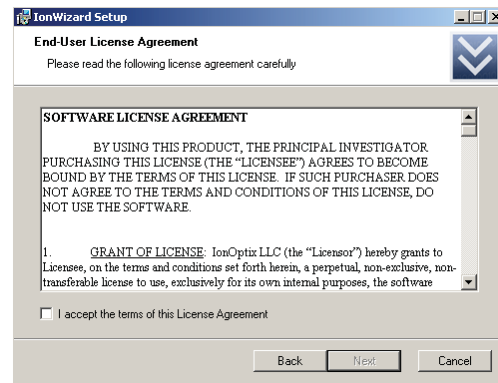


If this is the first time that you have installed this version of IonWizard you will see this dialog.

Click the Migrate Current button to copy the configuration files from the current installation.

Click the Use Blank button to set the configuration for this version to blank (empty) configuration files.



If you chose Use Blank the configuration files from the current version will NOT be erased - they will just not be copied to the new version.

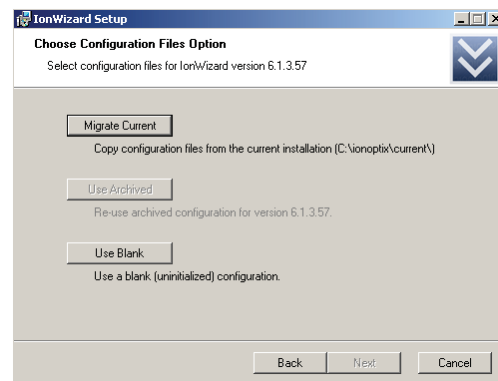

If you have previously installed a version of IonWizard, removed it and are now installing the same version again you will see this dialog.

Click the Migrate Current button to copy the configuration files from the current installation.

Click the Use Archived button to use the configuration files that that were last used with this version of IonWizard

Click the Use Blank button to overwrite the archived configuration files for this version with blank (empty) configuration files.



Migrate Current and Use Blank will replace the archived configuration for the version that is being installed. Configuration files for other versions will NOT be affected.




This screen is you final screen that you see before upgrade begins. If you need to review your choices you can **click** the Back button to return to previous screens or **click** the Cancel button to exit the installer.

Click the Upgrade button when you are ready to start the upgrade


As IonWizard is upgarding you will see this progress screen. The Status text ("Copying new files" in the sample picture) will change to indicate the current operation and the bar underneath will animate to show progress.



The progress bar will fill multiple times during installation.

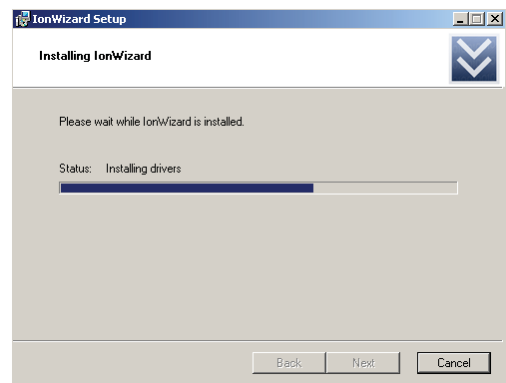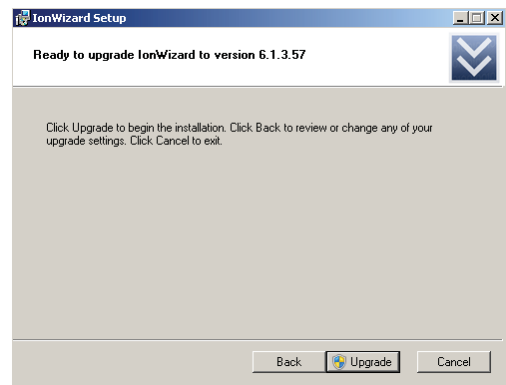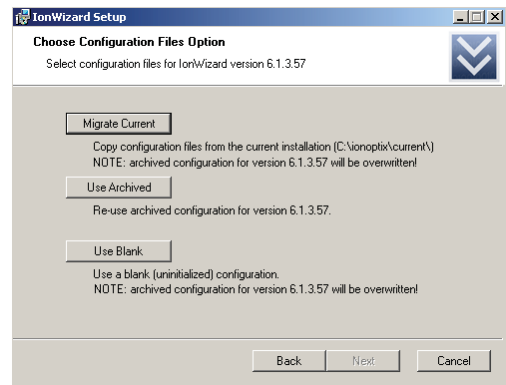

If the installation was completed without error then you will see this dialog. **Click** the **Finish** *button* to exit the installer. You may then run IonWizard using the Windows Start Menu. See [Start menu and Control Panel options](#) ¹³

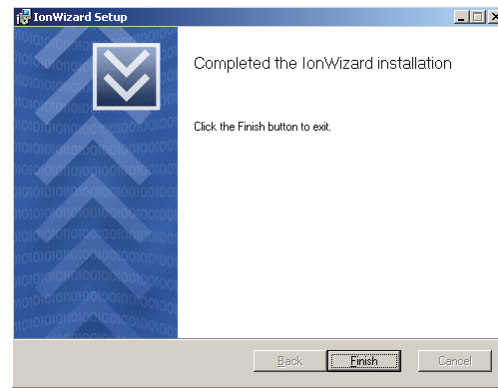

If an error is encountered during upgrade the installer will stop and display this message. **Click** the **Finish** *button* to exit the installer.

If you encounter an error please try running the installer a second time **AFTER** restarting Windows before contacting support for help.

